MANUAL


Lab Vac LVH40

(GB) Installation
(ES) Instalación
(SE) Installation

(DK) Installation
(FR) Installation
(FI) Asennus

(DE) Einbau
(IT) Installazione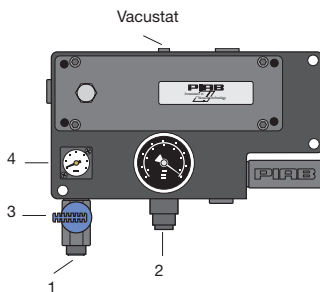

(GB)

1. Connect compressed air to port 1.
2. Connect vacuum hose (not included in delivery).
3. Start the pump.
4. Adjust pressure if necessary (0.4-0.6 MPa, 58-87 psi).

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Compressed air recommendations

Piab recommends filtered, oil-free compressed air for its multi-stage ejectors.

Maximum particle size:	40 μ
Maximum particle concentration:	10 mg/m ³ (atmospheric air)
Maximum oil content:	1 mg/m ³ (atmospheric air)

Compressed air supply

Piab LVH40 is designed for a compressed air supply pressure of 0.4-0.6 MPa (58-87 psi). To ensure that the supply pressure to the pump will be correct, it is important for the compressed air lines to be correctly sized. Use at least the sizes specified in the table below.

Hose length 1 m	Hose length 2 m	Hose length 3 m
Ø 5.2 mm	Ø 6.0 mm	Ø 8.2 mm
<i>The dimensions are calculated for a pressure drop of 0.01 MPa (0.1 bar).</i>		

Insufficient feed pressure

Piab LVH40 will not be damaged in case of insufficient feed pressure. What happens is that the flow decreases, and the maximum vacuum level decreases according to the table below.

Feed pressure, MPa	Max. vacuum level, mbar abs
0.60	20
0.55	20
0.50	20
0.45	50
0.40	270
0.35	420
0.30	550

Resistance to chemicals

Chemicals	Material in pump	Recommended seal and flap valve material		
		EPDM seal	Viton® (FPM) seal	Kalrez®
Acetone nitrile	A	A	U	A
Ammonia	A	A	U	A
Hydrofluoric acid	A	U	A	A
NaOH	A	A	B	A
Nitric acid	B	A	A	A
Hydrochloric acid	A	A	A	A
Sulphuric acid	A	B	A	A

*PPS = Polyphenylene sulphide
A = Recommended – little or no effect
B = Minor to moderate effect
C = Moderate to severe effect
U = Not recommended

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(GB)

- Check the feed pressure and blockage or leakage in the vacuum hose.
- Disconnect compressed air and dismantle pump house. Take care to check the sequence of assembly and position of gaskets and flap valves.
- Clean pump house. Check and clean filter, gaskets and flap valves thoroughly.
- Replace parts in reversed order. Note! The position of flap valves.

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